


Acetic Acid monitoring in Ethanol



Column: **MEGA-FFAP EXT FAST - 0.15mm, 0.25 μ m, 10m**

Catalog Code: F-FFAPEXT-015-025-10

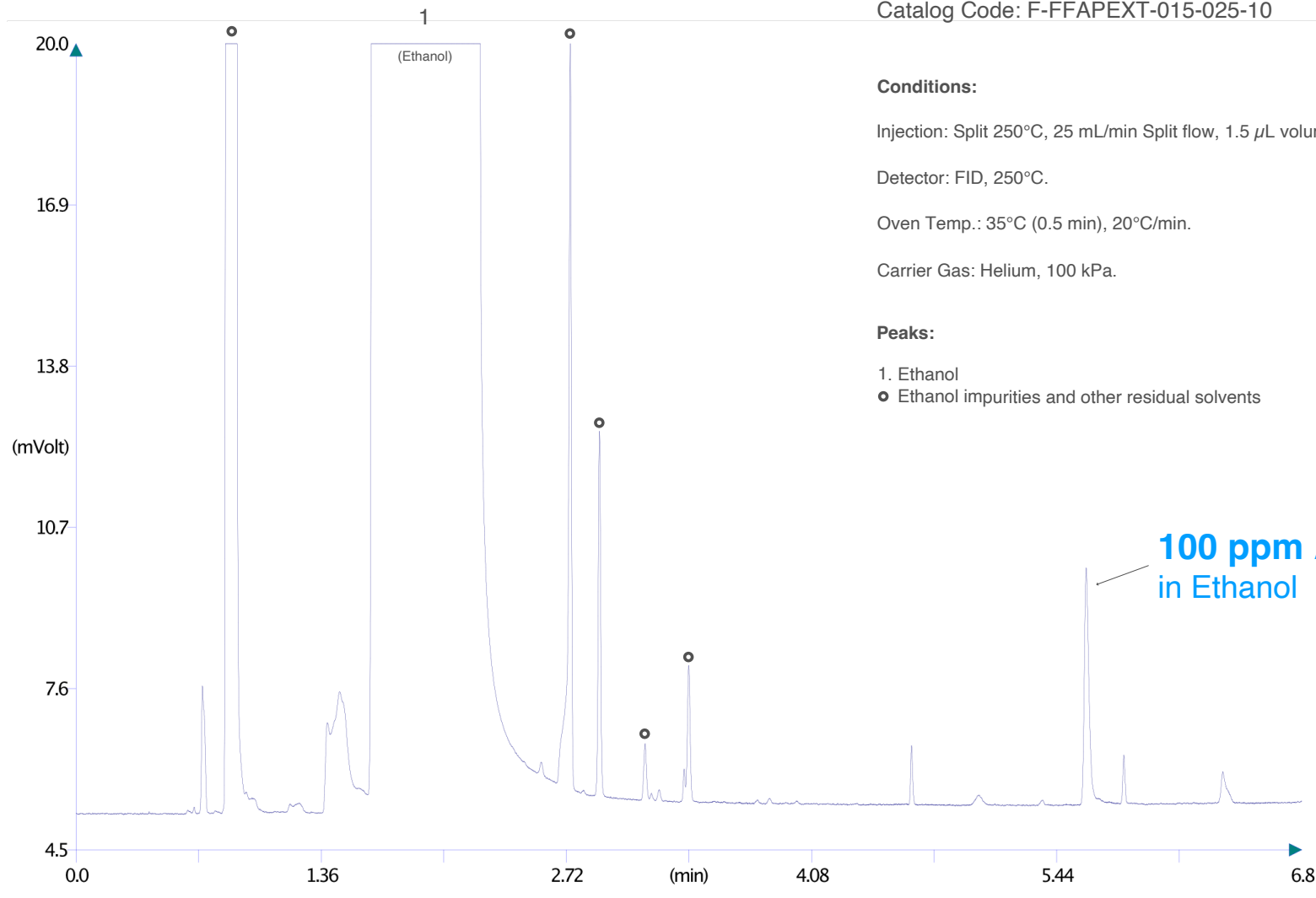

Conditions:

Injection: Split 250°C, 25 mL/min Split flow, 1.5 μ L volume injected.

Detector: FID, 250°C.

Oven Temp.: 35°C (0.5 min), 20°C/min.

Carrier Gas: Helium, 100 kPa.

Peaks:

- 1. Ethanol
- Ethanol impurities and other residual solvents