


Column: **MEGA-1701 - 0.32mm, 1.00 μ m, 30m**

Catalog Code: S-1701-032-100-30

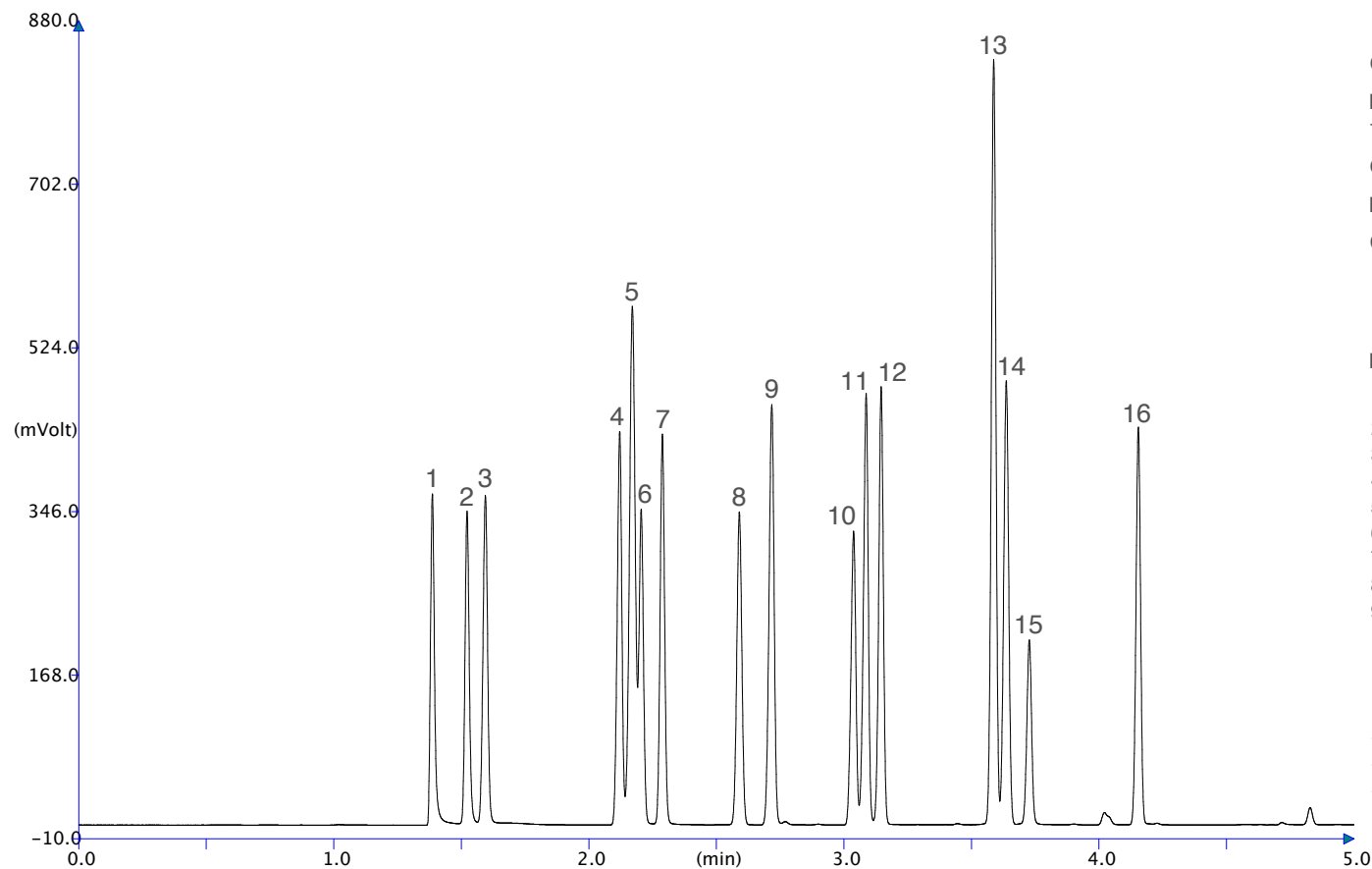

Conditions:

Injection: Split 250°C, 0,2 μ L (with micro-syringe), 150mL/min Split Flow.

Oven Temp.: 35°C (0.5 min), 20°C/ min, 250°C.

Detector: FID, 280°C.

Carrier Gas: Hydrogen, 60 kPa.

Peaks:

1. Ethanol
2. Acetone
3. 2-propanol
4. 1-propanol
5. Isobutanol
6. Ethyl acetate
7. Methyl-Ethyl-Ketone
8. Isopropyl acetate
9. Cyclohexane
10. Methoxy propanol
11. 1-butanol
12. Propyl acetate
13. Toluene
14. Isobutyl-Methyl-Ketone
15. Ethoxy propanol
16. Butyl acetate