

Nerolidol (+/-, *cis*, *trans*)


Chiral separation on MEGA-DEX columns

CHIRAL DEX

MEGA[®]
gas chromatography solutions

Column: **MEGA-DEX DET Beta - 0.25mm, 0.25 μ m, 25m**

Catalog Code: DEX-DET-B-025-025-25

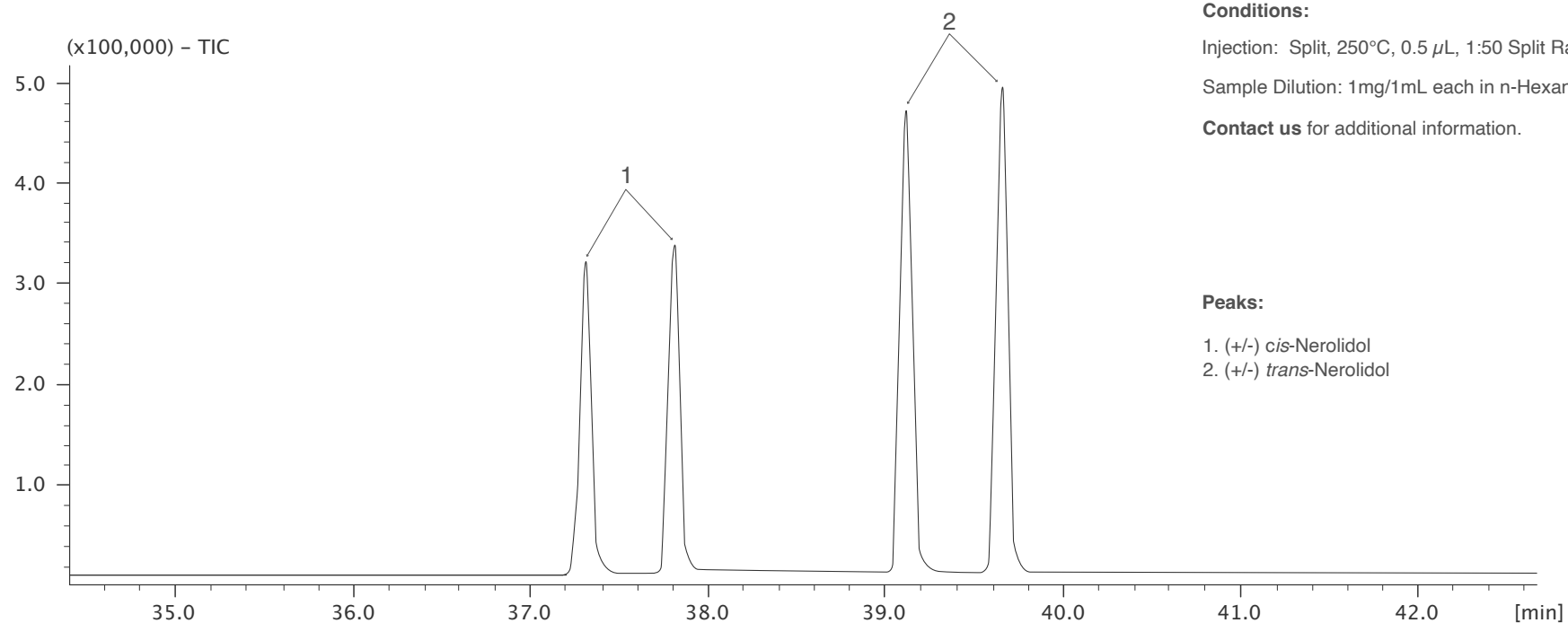

Conditions:

Injection: Split, 250°C, 0.5 μ L, 1:50 Split Ratio.

Sample Dilution: 1mg/1mL each in n-Hexane.

Contact us for additional information.

Peaks:

1. (+/-) *cis*-Nerolidol
2. (+/-) *trans*-Nerolidol

Acknowledgement: Prof C. Bicchi, C. Cordero et al. - Università di Torino, Dipartimento di Scienza e Tecnologia del Farmaco, Via Pietro Giuria, 9 - Torino - Italy.