

Alkanes C₈ - C₁₂ + Aromatics C₆ - C₁₀

CUSTOM
DEDICATED

MEGA[®]
gas chromatography solutions

Column: **MEGA-TCEP - 0.25mm, 0.40μm, 60m**

Catalog Code: C-TCEP-025-040-60

Conditions:

Injection: Split, 200°C.

Detector: FID, 200°C.

Oven Temp.: 95°C (5 min), 2.5°C/min.

Carrier Gas: Nitrogen, 90 kPa.

Peaks:

1. n-octane
2. n-nonane
3. n-decane
4. n-undecane
5. n-dodecane
6. Benzene
7. Toluene
8. Ethylbenzene
9. m-xylene
10. p-xylene
11. n-propylbenzene
12. o-xylene
13. n-butylbenzene

